



 Product Information
 ExcelBand™ series

 3-color Broad Range Protein Marker,
 (3.5-245 kDa)

## PM2700

Size:  $250 \ \mu l \times 2$ 5 \ \mu l per well for 100 applications

#### Storage

4°C for 3 months -20°C for 24 months

#### Loading Volume

- 3 µl or 5 µl per loading for clear visualization during electrophoresis on a 15-well or 10-well mini-gel.
- **1.5~2.5 µl** per well for general Western blotting transfer.
- Apply more for a thicker (> 1.5 mm) or larger gel.

## Description

The PM2700 ExcelBand<sup>™</sup> 3-color Broad Range Protein Marker is a ready-to-use three-color protein standard with 13 pre-stained proteins covering a wide range of molecular weights from 5 to 245 kDa in Tris-Glycine buffer (3.5 kDa to 235 kDa in Bis-Tris (MOPS) buffer and Bis-Tris (MES) buffer). Proteins are covalently coupled with a blue chromophore except for two reference bands (one green and one red band at 25 kDa and 75 kDa, respectively) when separated on SDS-PAGE (Tris-Glycine buffer).

The PM2700 ExcelBand<sup>™</sup> 3-color Broad Range Protein Marker is designed for monitoring protein separation during SDS-polyacrylamide gel electrophoresis, verification of Western transfer efficiency on membranes (nitrocellulose, PVDF, or nylon) and for approximating the size of proteins.

The marker is supplied in gel loading buffer and is ready to use. Do NOT heat, dilute, or add reducing agents before loading.

## Contents

Approximately 0.1 $^{0.4}$  mg/ml of each protein in the buffer (20 mM Tris-phosphate (pH 7.5 at 25 $^{\circ}$ C), 2 % SDS, 0.2 mM Dithiothreitol, 3.6 M Urea, and 15 % (v/v) Glycerol).

## **Guide for Molecular Weight Estimation**

Migration patterns and approximate MWs (kDa) of PM2700 ExcelBand<sup>™</sup> 3-color Broad Range Protein Marker in different electrophoresis conditions are listed below:

Band	Color	TRIS- GLYCINE	BIS-TRIS (MOPS)	BIS-TRIS (MES)	
1	Blue	245	235	235	
2	Blue	180	170	170	
3	Blue	140	130	130	
4	Blue	100	93	93	
5	Red	75	70	72	
6	Blue	60	53	53	
7	Blue	45	41	42	
8	Blue	35	30	30	
9	Green	25	22	23	
10	Blue	20	18	18	
11	Blue	15	14	14	
12	Blue	10	9	10	
13	Blue	~5	~3.5	~3.5	



# ExcelBand™ 3-color Broad Range Protein Marker PM2700



Note. The apparent molecular weight (kDa) of each protein has been determined by calibration against an unstained protein standard; supplemental data should be considered for more accurate adjustment in different electrophoresis conditions.

## **Quality Control**

Under suggested conditions, PM2700 ExcelBand<sup>™</sup> 3color Broad Range Protein Marker resolves 13 major bands in SDS-PAGE (Tris-Glycine, MOPS, and MES buffer) and after Western blotting transfer to a nitrocellulose membrane.

## Other information

SMOBIO Technology, Inc. claims all warranties with respect to this document, expressed or implied, including but not limited to those of merchantability or fitness for a particular purpose. In no event shall SMOBIO Technology, Inc. be liable, whether in contract, tort, warranty, or under any statute or any other basis for special, incidental, indirect, punitive, multiple or consequential damages in connection with or arising from this document, including but not limited to the use thereof.

Caution: Not intended for human or animal diagnostic or therapeutic uses.