
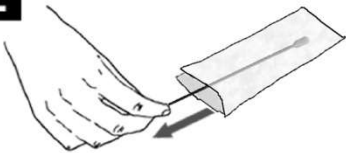





## Instructions for use of Isohelix RD-01 Buccal Swabs with RapiDri Pouch

REF RD-01

Intended for the retrieval of buccal cells. Single use only  
Store at room temperature



<p>Take your DNA sample at least one hour after eating, drinking or cleaning your teeth. For best results, rinse mouth with water immediately prior to sampling.</p>	
<p><b>1</b></p> 	<p>Pull open the package from one end.</p>
<p><b>2</b></p> 	<p>Remove the swab from the wrapper, taking care not to touch the white swab head with your fingers.</p>
<p><b>3</b></p> 	<p>Insert the swab into your mouth and rub firmly against the inside of your cheek or underneath lower or upper lip. For standard DNA collection rub for <b>1 minute</b> and in all cases rub for a <b>minimum</b> of 20 seconds. <b>Important – use reasonable, firm and solid pressure</b></p>
<p><b>4</b></p> 	<p>Place the swab head first into the RapiDri pouch so that the head sits within the drying area.</p>
<p><b>5</b></p> 	<p>Fold the RapiDri Pouch at the Fold Line, with the clear side inwards as shown, and use the barcoded sticker to seal.</p>