# **Anti-Conjugated Taurine**

# Product Information Sheet # 1021GE



#### **SUMMARY**

shipped on blue ice; store at -20 °C

For research use only.

Not for use in diagnostic or therapeutic procedures.

### **Description**

Polyclonal antisera were raised in rabbits after immunization with the conjugates: Taurine-glutaraldehyde-carriers.

**Contents:** The antibody is supplied in a unit size of 50 µl in PBS with 50% glycerol.

## Storage and Handling

When this product is stored undiluted at -20 °C or lower, it is stable for at least 2 years. Multiple freezing and thawing is only possible up to 5 times. The antibodies can be prediluted 10x in PBS containing 0.1% merthiolate or a mixture of PBS/glycerol (1:1, vol./vol.). At 4 °C these solutions are stable for 2 months.

## Specificity

Using a conjugate Taurine-glutaraldehyde-protein, antibody specificity was determined in an ELISA test by competition experiments with the following compounds:

Compound	Cross-	Cross-reactivity ratio (a)	
Taurine-G-PL	1		
Hypotaurine-G-PL	1:	15	
GABA-G-PL	1:	560	
p-Alanine-G-PL	1:	1,860	
Aspartate-G-PL	1:	5,040	
Taurine=G=PL (b)	1:	6,700	
Glycine-G-PL	1:	14,200	
Cysteine-G-PL	1:	> 20,000	
Glutamate-G-PL	1:	> 20,000	
Taurine	1:	> 40,000	

(a): Taurine-G-PL concentration/unconjugated or conjugated aminoacids concentration at half displacement;

#### **Recommended dilution**

The anti-conjugated Taurine antibodies can be diluted 1:2,000 to 1:5,000.

MoBiTec GmbH, Germany Phone: +49 551 70722 0 Fax: +49 551 70722 22 E-Mail: info@mobitec.com www.mobitec.com

<sup>(</sup>b): Non-reduced conjugate

G = Glutaraldehyde, PL = Poly-lysine

# **Anti-Conjugated Taurine**

# Product Information Sheet # 1021GE



#### Reference

CAMPISTRON G., GEFFARD M. and BUIJS R.M., Immunological approach to the detection of taurine and immunocytochemical results. J. Neurochem. (1986) 46, 862-868.

# **Order Information, Shipping and Storage**

Order#	Product	Quantity
1021GE	Anti-Conjugated Taurine	50 μl
shipped on blue ice; store at -20 °C		

## **Contact and Support**

MoBiTec GmbH ● Lotzestrasse 22a ● D-37083 Goettingen ● Germany

Customer Service – General inquiries & orders Technical Service – Product information

phone: +49 (0)551 707 22 0 phone: +49 (0)551 707 22 70 fax: +49 (0)551 707 22 22 fax: +49 (0)551 707 22 77 e-mail: order@mobitec.com e-mail: info@mobitec.com

MoBiTec in your area: Find your local distributor at www.mobitec.com

MoBiTec GmbH, Germany Phone: +49 551 70722 0 Fax: +49 551 70722 22 E-Mail: info@mobitec.com www.mobitec.com