# **Anti-Conjugated Spermine**

# Product Information Sheet # 1047GE



### **SUMMARY**

shipped on blue ice; store at -20 °C

For research use only.

Not for use in diagnostic or therapeutic procedures.

# **Description**

Polyclonal antisera were raised in rabbits after immunization with the conjugates: Spermineglutaric anhydride-carriers.

**Contents:** The antibody is supplied in a unit size of 50 µl in PBS with 50% glycerol.

### Storage and Handling

When this product is stored undiluted at -20 °C or lower, it is stable for at least 2 years. Multiple freezing and thawing is only possible up to 5 times. The antibodies can be prediluted 10x in PBS containing 0.1% merthiolate or a mixture of PBS/glycerol (1:1, vol./vol.). At 4 °C these solutions are stable for 2 months.

## **Specificity**

Using a conjugate Spermine-Glutaric Anhydride-protein, antibody specificity was determined in an ELISA test by competition experiments with the following compounds:

Compound	Cross-reactivity ratio (a)	
Spermine-GA-BSA	1	
Spermidine-GA-BSA	1:3	
Diamino Pentane-G-BSA	1:>700	
Ornithine-GA-BSA	1:>500	

<sup>(</sup>a): Spermine-GA-BSA concentration / concentration of unconjugated or conjugated polyamine at half displacement;

#### **Recommended dilution**

The anti-conjugated Spermine antibodies can be diluted 1:2,000 to 1:5,000.

MoBiTec GmbH, Germany Phone: +49 551 70722 0 Fax: +49 551 70722 22 E-Mail: info@mobitec.com www.mobitec.com

G = Glutaraldehyde, GA = Glutaric Anhydride, BSA = Bovine Serum Albumine

# **Anti-Conjugated Spermine**

# Product Information Sheet # 1047GE



# **Order Information, Shipping and Storage**

Order#	Product	Quantity
1047GE	Anti-Conjugated Spermine	50 µl
shipped on blue ice; store at -20 °C		

## **Contact and Support**

MoBiTec GmbH ● Lotzestrasse 22a ● D-37083 Goettingen ● Germany

Customer Service – General inquiries & orders Technical Service – Product information

phone: +49 (0)551 707 22 0 phone: +49 (0)551 707 22 70 fax: +49 (0)551 707 22 22 fax: +49 (0)551 707 22 77 e-mail: order@mobitec.com e-mail: info@mobitec.com

MoBiTec in your area: Find your local distributor at www.mobitec.com

MoBiTec GmbH, Germany Phone: +49 551 70722 0 Fax: +49 551 70722 22 E-Mail: info@mobitec.com www.mobitec.com