

AquaGenomic Instruction Manual

General Information

Description

AquaGenomic™ is a multifunctional aqueous reagent for genomic DNA extraction. It may be used to extract DNA from all types of specimens, from bacteria to plant and animal tissues. The protocols are simple, fast, and scalable. AquaGenomic is nontoxic and its lysate may be used for PCR directly without DNA purification. AquaGenomic can efficiently extract DNA from dried specimen swabs. It may facilitate the use of dried specimen swabs for low-cost biobanking, biosurveillance and epidemiological studies.

Specification

Product Name	AquaGenomic™ Solution
Product #	2001, 2030
Size	2001: 1 ml AquaGenomic for 10 minipreps from cultured cells 2030: 30 ml AquaGenomic for 300 minipreps from cultured cells
MSDS	Available at www.multitargetpharm.com
Storage	Store tightly capped at room temperature (~22°C).

Terms & Condition

Product Usage: For In Vitro Laboratory Research Use Only. NOT to be administered to humans or used for medical diagnosis.

Warranties and Liabilities: MultiTarget Pharmaceuticals accepts no responsibility and shall not be held liable for any loss, damage, expense, consequential, or accidental damage, including damage to property, person, or premises arising out of the use, the results of use, or the inability to use these products. MultiTarget Pharmaceuticals MAKES NO WARRANTIES, EXPRESSED OR IMPLIED, INCLUDING, WITHOUT LIMITATION, WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR USE.

Patents, Trademarks & Copyrights

AquaGenomic™ is a trademark of MultiTarget Pharmaceuticals, LLC.
©2018 Multitarget Pharmaceuticals, LLC. All rights reserved.

AquaGenomic Cell Protocol

This protocol can be used to prepare 5-10 µg of genomic DNA from 1-2 million cultured cells. For other sample sizes, use additional 100 µl of AquaGenomic Solution for each million nucleated cells.

1. Harvest the Cells

Pellet ~0.5-2 million cultured cells in a 1.5-ml microfuge tube by centrifugation at 12,000 xg for 60 sec. Aspirate or decant to discard the supernatant. (*Note: For convenience, cell culture may be mixed with 1 vol of AquaGenomic for DNA extraction without needing to pellet the cells.*)

2. Extract the DNA

Add 100 µl of AquaGenomic to the cell pellet. Suspend and lyse the cells by vortex vigorously for 30-60 sec. Incubate at room temperature or at 75° C for 20 min (75° C incubation produces better lysis and inactivates any residual DNases). (*Note: To use the lysate for PCR, dilute an aliquot of the lysate with 10-20 vol of water and then use 0.25 µl of the diluted lysate in a 25-µl PCR reaction.*)

3. Pellet the DNA

Add 1 vol of isopropanol (e.g., to 100 µl crude lysate, add 100 µl isopropanol) and vortex for 30-60 sec to mix well. Centrifuge at 12,000 xg for 5 min to pellet the DNA. Decant to discard the supernatant.

4. Rinse the DNA

Gently fill up the tube and its lid with 70% ethanol from a squirt bottle and then flip the tube to discard the ethanol solution. Repeat the 70% ethanol rinse once. Flip the tube forcefully a few times and tap it on a paper towel to remove residual ethanol. Let the DNA pellet air-dry for 5-10 min.

5. Solubilize the DNA

Add 100 µl of TE buffer or deionized water to the DNA pellet, pipette or vortex vigorously to suspend the DNA. Incubate at 22 °C for 15 min (*Optional: To fully solubilize and recover the DNA, incubate at 65-75 °C for 15 min or at room temperature overnight*). Centrifuge at 12,000 xg for 5 min to pellet the insoluble and transfer the clear DNA solution to a new tube.

AquaGenomic Tissue Protocol

This protocol may be used to extract DNA from tissues (including common research specimens, such as drosophila, mouse-tail snip, nematode, zebrafish, and plant tissues) by homogenization or by Proteinase K digestion in AquaGenomic solution.

1. Harvest the Cells

Cut out a ~2 mm cube (~10 mg) of frozen or fresh tissue.

2. Extract the DNA

By homogenization: Homogenize the tissue in 200 µl of AquaGenomic. Move the pestle up and down slowly while vortexing at top speed to enhance homogenization. Alternatively, homogenize the tissue in 0.5-ml screw-capped tubes with a multichannel bead beater. Transfer the homogenate to a 1.5-ml microfuge tube. Incubate at room temperature or at 75° C for 20 min (75° C incubation produces better lysis and inactivates any residual DNases).

By Proteinase K digestion: Place the tissue (e.g., a mouse tail clip) in a microfuge tube preloaded with 200 µl of AquaGenomic containing 10 µg of Proteinase K. Incubate at 55 °C for >90 min to digest the tissue, and then at 95 °C for 10 min to inactivate the Proteinase K. The tissue is readily disintegrated by vortexing or pipetting.

3. Pellet the debris

Centrifuge at 12,000 xg for 5 min to pellet the debris. (*Note: To use the lysate for PCR, dilute an aliquot of the lysate with 10-20 vol of water and then use 0.25 µl of the diluted lysate in a 25-µl PCR reaction.*)

4. Pellet the DNA

Transfer the clear lysate (~180 µl) to a new 0.5-ml microfuge tube. Add 1 vol (~180 µl) of isopropanol and vortex for 60 sec to mix well. Centrifuge at 12,000 xg for 5 min to pellet the DNA. Decant to discard the supernatant. Gently fill up the tube and its lid with 70% ethanol from a squirt bottle and then flip the tube to discard the ethanol solution. Repeat the 70% ethanol rinse once. Flip the tube forcefully a few times and tap it on a paper towel to remove residual ethanol. Let the DNA pellet air-dry for 5-10 min. Add 100 µl of TE buffer or deionized water to the DNA pellet, pipette or vortex vigorously to suspend the DNA. Incubate at 22 °C for 15 min (*Optional: To fully solubilize and recover the DNA, incubate at 65-75 °C for 15 min or at room temperature overnight*). Centrifuge at 12,000 xg for 5 min to pellet any insoluble and transfer the DNA solution to a new tube.

AquaGenomic Saliva Protocol

Approximately 10-20 µg of genomic DNA can be obtained from 100 µl of saliva or a buccal swab or 200 µl of mouthwash, using one of the following methods.

1. Harvest the Cells

- (a) Saliva. Swirl and rub your tongue against the inside of your cheek and gum for ~5-10 times. Carefully spit the saliva into a clean weight boat or a 15-ml conical tube.
- (b) Swab. Use a swab to rub the inside of your cheek and gum for ~5-10 times and let it soak up the saliva. Air-dry the swab in its pouch.
- (c) Mouthwash. Swirl and rub your tongue against the inside of your cheek and gum for ~5-10 times. Rinse the mouth with 10-20 ml of Scope mouthwash and spit it into a 50-ml conical tube.

2. Extract the DNA

- (a) Saliva. Transfer 100 µl of saliva to a microfuge tube preloaded with 100 µl of AquaGenomic solution and incubate at room temperature or at 75° C for 20 min (75° C incubation produces better lysis and inactivates any residual DNases).
- (b) Swab. Cut off the tip of the swab into a 1.5-ml microfuge tube. Add 300 µl of AquaGenomic solution and incubate at 75 °C for 20 min. Use a 1-ml pipet tip to smash the swab 10 times to the bottom of the tube to squeeze out of the liquid.
- (c) Mouthwash. Centrifuge 200 µl of mouthwash at 10,000 xg for 5 min to pellet the buccal cells and discard the supernatant. Add 200 µl of AquaGenomic solution. Vortex to mix well and incubate at 22 °C for 10 min.

(Note: To use the lysate for PCR, dilute an aliquot of the lysate with 10-20 vol of water and then use 0.25 µl of the diluted lysate in a 25-µl PCR reaction.)

3. Pellet the DNA

Add 1 vol of isopropanol to the crude lysate containing the cell debris (e.g., to 200 µl crude lysate, add 200 µl isopropanol) and vortex for 60 sec to mix well. Centrifuge at 12,000 xg for 5 min to pellet the DNA. Decant to discard the supernatant. Gently fill up the tube and its lid with 70% ethanol from a squirt bottle and then flip the tube to discard the ethanol solution. Repeat the 70% ethanol rinse once. Flip the tube forcefully a few times and tap it on a paper towel to remove residual ethanol. Let the DNA pellet air-dry for 5-10 min. Add 100 µl of TE buffer or deionized water to the DNA pellet, pipette or vortex vigorously to suspend the DNA. Incubate at 22 °C for 15 min (*Optional: To fully solubilize and recover the DNA, incubate at 65-75 °C for 15 min or at room temperature overnight*). Centrifuge at 12,000 xg for 5 min to pellet the insoluble, and transfer the clear DNA solution to a new tube.

AquaGenomic Microbe Protocol

This protocol can be used to prepare 10-20 µg of DNA from 1 ml overnight microbial culture. For other preparation scales, use 200 µl of AquaGenomic Solution for each milliliter of overnight culture.

1. Harvest the Cells

Centrifuge 1 ml overnight bacterial culture at 12,000xg for 60 sec to pellet the cells. Aspirate or decant to discard the supernatant.

2. Extract the DNA

For Gram-negative bacteria: Add 200 µl of AquaGenomic Solution to the cell pellet. Suspend the cells by vortexing vigorously for 30 sec. Alternatively, homogenize the samples in AquaGenomic in 0.5-ml screw capped tubes with a multichannel bead beater.

For Gram-positive bacteria or yeast: Treat the bacterial or yeast cells with lysozyme or lyticase (not supplied) according the enzyme manufactures' instruction. Add ~20 mg of sand (Sigma # 274739, white, 50+70 mesh) and 200 µl of AquaGenomic Solution containing 100 µg/ml Proteinase K to the sample. Incubate at 55° C for 60 min and then at 95 °C for 10 min to inactivate the Proteinase K.

(Note: To use the lysate for PCR, dilute an aliquot of the lysate with 10-20 vol of water and then use 0.25 µl of the diluted lysate in a 25-µl PCR reaction.)

3. Pellet the DNA

Add 1 vol of isopropanol to the crude lysate containing the cell debris (e.g., to 200 µl crude lysate, add 200 µl isopropanol) and vortex for 30-60 sec to mix well. Centrifuge at 12,000 xg for 5 min to pellet the DNA and cell debris. Decant to discard the supernatant. Gently fill up the tube and its lid with 70% ethanol from a squirt bottle and then flip the tube to discard the ethanol solution. Repeat the 70% ethanol rinse once. Flip the tube forcefully a few times and tap it on a paper towel to remove residual ethanol. Let the DNA pellet air-dry for 5-10 min. Add 100 µl of TE buffer or deionized water to the DNA pellet, pipette or vortex vigorously to suspend the DNA. Incubate at 22 °C for 15 min (*Optional: To fully solubilize and recover the DNA, incubate at 65-75 °C for 15 min or at room temperature overnight*). Centrifuge at 12,000 xg for 5 min to pellet the insoluble, and transfer the clear DNA solution to a new tube.

AquaGenomic Stool and Soil Protocol

This protocol uses 200 µl of AquaGenomic Solution to prepare 5-10 µg of DNA from 15 mg of feces. AquaPrecipi solution (Item # 3015, not included) is generally required to purify fecal and soil DNA and remove PCR inhibitors. However, for mouse fecal DNA extraction and genotyping, AquaPrecipi purification may not be necessary.

1. Harvest the Cells

Weigh out 15 mg of wet feces (~10 mg of dry fecal pellet or a mouse fecal pellet) or 30 mg of soil in a 1.5-ml microfuge tube.

2. Extract the DNA

Add 200 µl of AquaGenomic solution to the sample. Incubate at room temperature or at 75 °C for 20 min (75 °C incubation produces better lysis and inactivates any residual DNases). Homogenize the sample with a microfuge pestle or vortex vigorously for 1-2 min. Alternatively, homogenize the samples in AquaGenomic in 0.5-ml screw capped tubes with a multichannel bead beater. If mitochondrial DNA extraction is desired, add Proteinase K to AquaGenomic to 100 µg/ml, incubate at 55 °C for 60 min to digest the mitochondria and then at 95 °C for 15 min to inactivate the Proteinase K.

3. Pellet the Debris

Vortex vigorously for 60 sec and centrifuge at 12,000 xg for 5 min to pellet the debris. Transfer the clear lysate (~100 µl) to a new 0.5-ml microfuge tube (*Note: The lysates of most fecal and soil samples cannot be used directly in PCR reactions as they contain large amounts of PCR inhibitors. AquaPrecipi is required for the removal of these fecal and soil PCR inhibitors in the next step.*).

4. Pellet the DNA

Add 0.5 vol (~50 µl) of AquaPrecipi (#3015, order separately) and 0.5 vol (~50 µl) of ethanol. Vortex for 60 sec and centrifuge at 12,000 xg for 5 min to pellet the DNA. Decant to discard the supernatant. Gently fill up the tube and its lid with 70% ethanol from a squirt bottle and then flip the tube to discard the ethanol solution. Repeat the 70% ethanol rinse once. Flip the tube forcefully a few times and tap it on a paper towel to remove residual ethanol. Let the DNA pellet air-dry for 5-10 min. Add 100 µl of TE buffer or deionized water to the DNA pellet, pipette or vortex vigorously to suspend the DNA. Incubate at 22 °C for 15 min (*Optional: To fully solubilize and recover the DNA, incubate at 65-75 °C for 15 min or at room temperature overnight*). Centrifuge at 12,000 xg for 5 min to pellet any insoluble, which contains residual PCR inhibitors, and transfer the clear DNA solution to a new tube.

AquaGenomic Swab Protocol

Cotton swabs are commonly used in forensic evidence collection. However, existing DNA extraction methods can recover only ~100-500 ng of DNA from a dried specimen swab as the majority of DNA remains entrapped within the cotton matrix. With AquaGenomic, ~5 µg of DNA can be extracted from a dried specimen swab. It is possible to use cotton swabs for specimen collection, transportation, storage, and DNA extraction for low-cost biobanking, biosurveillance, and epidemiological studies.

1. Collect the specimen. Use a cotton swab to soak up the specimen (~100-200 µl), such as blood, saliva, mucus, semen, feces, cultured mammalian or bacteria cells, homogenized animal or plant tissues, or any other potential sources of biospecimens. Air-dry the specimen swab at ambient temperature (20-50 °C) for >24 hours. The dried specimen swabs can be shipped at ambient temperature and then stored at room temperature in sealed paper envelopes or plastic bags with desiccant for many years.

2. Extract the proteins and small molecules (dried blood swabs). Cut off the specimen swab tip into a 1.5-ml microfuge tube. Add 400 µl deionized water or buffer. Soak the swab for >30 min. Use a 1-ml pipet tip to smash the swab ~10 times and press it to the bottom of the tube to squeeze out the solution. Transfer as much liquid as possible (~300 µl) to a new 1.5-ml microfuge tube for analysis of proteins or small molecules.

3. Extract the DNA. Add 300 µl AquaGenomic to the 1.5-ml tube containing the wet swab from Step 2. Incubate at 22 °C for 1 hr and then at 75-85 °C for 20 min. Use a 1-ml blue pipet tip to smash the swab ~10 times and press it to the bottom of the tube to squeeze out the solution. Alternatively, homogenize the samples in AquaGenomic in 0.5-ml screw capped tubes with a multichannel bead beater. Transfer as much liquid as possible (~200 µl) to a new 0.5-ml microfuge tube or use a microfuge spin bucket to recover all the lysate from the swab to maximize DNA yield (*Note: To use the crude lysate for PCR, dilute an aliquot of the crude lysate with 10-20 vol of deionized water and use 0.25 µl of the diluted lysate in a 25-µl PCR reaction*).

4. Purify the DNA. Add 1 vol of isopropanol to the crude lysate (e.g., add 200 µl isopropanol to 200 µl lysate). Vortex to mix well and centrifuge at 12,000 xg for 5 min to pellet the DNA. Flip the tube to discard the isopropanol supernatant. Gently fill up the tube and its lid with 70% ethanol from a squirt bottle and then flip the tube to discard the ethanol solution. Repeat the 70% ethanol rinse once. Flip the tube forcefully a few times and tap it on a paper towel to remove residual ethanol. Let the DNA pellet air-dry for 5-10 min. Add 100 µl of TE buffer or deionized water to the DNA pellet, pipette or vortex vigorously to suspend the DNA. Incubate at 65-75 °C for 15 min or at room temperature overnight. Centrifuge at 12,000 xg for 5 min to pellet any insoluble and transfer the clear DNA solution to a new tube.

AquaGenomic Whole Blood Protocol

AquaGenomic may be used to extract total blood cellular and cell-free DNA from whole blood samples and its lysate may be used directly for PCR without further purification.

(a) To use lysate directly for PCR

1. Add 20 µl of AquaGenomic solution to each well in a 0.2-ml 96-well PCR plate.
2. Transfer 10 µl of well-mixed fresh or thawed whole blood to the AquaGenomic solution. Pipet up and down a few times to mix.
3. Incubate at 75 °C for 20 min in a PCR machine with heated lid. After the incubation, add 170 µl of water to the blood clot. Shake or pipet up and down a few times to mix.
4. To amplify the DNA, add 0.25 µl of the diluted lysate to 25 µl of PCR master mix and subject the reaction mix to 35-45 cycles of PCR amplification.

(b) To purify total blood DNA

1. Add 200 µl of AquaGenomic solution to a 0.5-ml tube.
2. Transfer 100 µl of well-mixed fresh or thawed whole blood to the AquaGenomic solution. Pipet up and down a few times to mix.
3. Incubate at 75 °C for 20 min. After the incubation, centrifuge the samples at 12,000 xg for 5 min. Transfer the clear lysate (~200 µl) to a new 0.5-ml microfuge tube.
4. Add 1 vol (~200 µl) of 100% isopropanol to the clear lysate and vortex for 60 sec to mix well. Centrifuge at 12,000 xg for 5-10 min at 22 °C to pellet the DNA.
5. Decant to discard the supernatant. Gently fill up the tube and its lid with 70% ethanol from a squirt bottle and then flip the tube to discard the ethanol solution. Repeat the 70% ethanol rinse once. Flip the tube forcefully a few times and tap it on a paper towel to remove residual ethanol. Let the DNA pellet air-dry for 5-10 min. Add 50 µl of TE buffer or deionized water to the DNA pellet, vortex vigorously to suspend the DNA (*Optional: To fully solubilize and recover the DNA, incubate at 65-75 °C for 15 min or at room temperature overnight*).

AquaGenomic WBC Protocol

This protocol can be used to prepare ~5-6 ug of genomic DNA from white blood cells (WBC) in 100 ul of whole blood. It requires a RBC Lysis Solution (Product # 4015, not included) to lyse and remove the RBC prior to DNA extraction from the WBC.

1. Harvest the Cells

Add 50 ul of RBC Lysis Solution to 100 ul fresh or thawed whole blood. Mix the contents by touch-vortex (a few seconds on and a split second off) for 20-30 times at top speed to lyse the RBC. Add 1 ml deionized water and vortex to wash the cells. Centrifuge at 12,000xg for 1 minute to pellet the WBC and nuclei. Aspirate to remove the supernatant.

2. Extract the DNA

Add 100 ul of AquaGenomic Solution to the cell pellet. Vortex to suspend the cell pellet. Incubate at room temperature or at 75° C for 20 min (75° C incubation produces better lysis and inactivates any residual DNases).

3. Pellet the Debris

Vortex vigorously for 30-60 seconds and centrifuge at 12,000xg for 5 minutes to pellet the debris. (*Optional: Add 5 ul isopropanol to the sample, vortex to reduce foaming and make the debris pellet become tighter following centrifugation*).

4. Pellet the DNA

Transfer the supernatant (~90 ul) to a new 0.5-ml microfuge tube. Add 1 vol (~90 µl) of isopropanol and vortex to mix well. Centrifuge at 12,000 xg for 5 min to pellet the DNA. Decant to discard the supernatant.

5. Rinse the DNA

Gently fill up the tube and its lid with 70% ethanol from a squirt bottle and then flip the tube to discard the ethanol solution. Repeat the 70% ethanol rinse once. Flip the tube forcefully a few times and tap it on a paper towel to remove residual ethanol. Let the DNA pellet air-dry for 5-10 min. Add 50 µl of TE buffer or deionized water to the DNA pellet, pipette or vortex vigorously to suspend the DNA. Incubate at 22 °C for 15 min (*Optional: To fully solubilize and recover the DNA, incubate at 65-75 °C for 15 min or at room temperature overnight*). Centrifuge at 12,000 xg for 5 min to pellet the insoluble and transfer the clear DNA solution to a new tube and store at -20 °C.

AquaGenomic Avian Blood Protocol

This protocol may be used to prepare about 50-100 µg of genomic DNA from 10 µl of avian blood sample stored in 100 µl of Queen's lysis buffer (10 mM Tris, 10 mM NaCl, 10 mM EDTA, 1% n-lauroylsarcosine, pH 7.5), using 200 µl of AquaGenomic solution and 100 µl of AquaRemove solution (order separately, #1208).

1. Collect the blood sample

Avian blood samples are routinely collected in Queen's lysis buffer at a ratio of 1 volume to 9-10 volumes, and stored at room temperature until DNA extraction.

2. Extract the DNA

Add 200 µl of AquaGenomic to 100 µl of lysed avian blood sample in Queen's lysis buffer. Vortex vigorously for 60 sec. Incubate at room temperature or at 75° C for 20 min (75° C incubation produces better lysis and inactivates any residual DNases).

3. Pellet the Debris

Add 100 µl of isopropanol diluted AquaRemove (order separately, #1208. *Note: Dilute the AquaRemove solution with an equal volume of isopropanol before use.*) and vortex vigorously for 60 sec to mix well. Centrifuge at 12,000 xg for 5 min to pellet the debris.

4. Pellet the DNA

Transfer the clear lysate (~400 µl) to a new 1.5-ml microfuge tube. Add 1 vol (~400 µl) of 100% isopropanol and vortex for 60 sec to mix well. Centrifuge at 12,000 xg for 5 min to pellet the DNA. Decant to discard the supernatant.

5. Rinse the DNA

Gently fill up the tube and its lid with 70% ethanol from a squirt bottle and then flip the tube to discard the ethanol solution. Repeat the 70% ethanol rinse once. Flip the tube forcefully a few times and tap it on a paper towel to remove residual ethanol. Let the DNA pellet air-dry for 5-10 min. Add 100 µl of TE buffer or deionized water to the DNA pellet, pipette or vortex vigorously to suspend the DNA. Incubate at 22 °C for 15 min (*Optional: To fully solubilize and recover the DNA, incubate at 65-75 °C for 15 min or at room temperature overnight*). Centrifuge at 12,000 xg for 5 min to pellet the insoluble and transfer the clear DNA solution to a new tube and store at -20 °C.

Frequently Asked Questions

Please read through these questions carefully. The answers provide additional helpful tips and useful information for the successful use of AquaGenomic.

1. Do I need to keep AquaGenomic in the freezer?

No, AquaGenomic Solution is stable at room temperature (~22 °C) for >1 year.

2. Does AquaGenomic Solution contain Proteinase K?

No. AquaGenomic can be used to extract DNA from most cells and tissues without needing protease digestion. However, adding Proteinase K (50 µg/ml) to AquaGenomic solution can increase DNA yield and is required for mitochondrial DNA extraction. You may homogenize the sample in AquaGenomic containing Proteinase K, incubate it at 55 °C for 1-2 hrs and then at 95 °C for 10-15 min to inactivate the Proteinase K.

3. I am worried about cross-contamination using homogenizers, any tips?

Between uses, you may wash the homogenizer with soap and running water, soak it in 10% bleach for ~5 min, and then rinse it with running deionized water. If you still feel uneasy, you may use Proteinase K digestion without using a homogenizer, or use a multichannel bead beater for homogenization in screw-capped tubes.

4. How may I maximize the DNA yield?

Incubating the DNA pellet in TE buffer or deionized water at 22 or 4 °C overnight before centrifuging to remove the insoluble may maximize DNA recovery.

5. Do I have to use the lysate immediately for PCR?

No, you may store the lysate at 4 °C until analysis. If the lysate has been incubated at 75-85 °C for 20 min, it may even be left at room temperature until analysis.

6. I got a weak PCR amplification using the lysate directly, how may I improve it?

You may try a few things to optimize the amplification: a) try to use different amount of lysate for the PCR, from 0.25 µl undiluted lysate to 20x diluted lysate, b) add 0.1 mg/ml BSA to the PCR reaction, c) add 1 mg/ml DTT to the PCR reaction, and d) increase the PCR cycle number to 45 cycles.

MATERIAL SAFETY DATA SHEET (MSDS)

TRADE NAME: AquaGenomic™ Solution
DATE OF ISSUE: March 1, 2015

SECTION I: PRODUCT AND MANUFACTURER INFORMATION

TRADE NAME: AquaGenomic™ Solution

MultiTarget Pharmaceuticals, LLC
5050 Edison Ave Ste 214, Colorado Springs, CO 80915, USA
Telephone: 1-801-769-6586
DATE PREPARED: 03-01-2015

SECTION II: COMPOSITION / INFORMATION ON INGREDIENTS

SYNONYMS: AquaGenomic™, genomic isolation solution
CHEMICAL CHARACTERIZATION: Proprietary aqueous solution of Sodium Dodocyl Sulfate and other non-hazardous chemicals.
HAZARDOUS COMPONENT: None.

SECTION III: HAZARD IDENTIFICATION

May be harmful if swallowed. May cause irritation. Avoid breathing vapors, or dusts. Avoid contact with eyes, skin, and clothes. Wash thoroughly after handling. Keep container closed.

SECTION IV: FIRST AID MEASURES

IF EYE CONTACT: Immediately flush eyes with copious amounts of water for at least 15 minutes. Assure adequate flushing of eyes by separating the eyelids with fingers.
IF SKIN CONTACT: Immediately wash skin with soap and copious amounts of water.
IF SWALLOWED: Wash out mouth with water provided person is conscious. Call a physician.

SECTION V: FIREFIGHTING MEASURES

EXTINGUISHING MEDIA: Use extinguishing media appropriate to surrounding fire conditions.
SPECIAL FIREFIGHTING PROCEDURES: Wear self-contained breathing apparatus and protective clothing to prevent contact with eyes and skin.

SECTION VI: ACCIDENTAL RELEASE MEASURES

PRECAUTIONARY MEASURES: Wear self-contained breathing apparatus, chemical safety goggles, rubber boots, and heavy rubber gloves.
CLEAN-UP PROCEDURES: Absorb on sand or vermiculite and place in closed container for disposal. Ventilate area and wash spill site after material pick-up is complete.

SECTION VII: HANDLING AND STORAGE

STORAGE: Store tightly closed at 20-25 °C.

SECTION VIII: EXPOSURE CONTROLS AND PERSONAL PROTECTION

Chemical safety goggles. Rubber gloves. Safety shower and eye bath. Wash thoroughly after handling. Do not get in eyes, on skin, or on clothing.

SECTION IX: PHYSICAL AND CHEMICAL PROPERTIES

PHYSICAL STATE / FORM: Liquid
COLOR: Colorless
ODOR: None
pH: (20 °C) N/V
VISCOSITY: (20 °C) N/A

MELTING POINT: N/A
BOILING POINT: N/A
IGNITION TEMPERATURE: N/A
FLASHPOINT: N/A
EXPLOSION LEVEL: N/A
VAPOR PRESSURE: (20 °C) N/A
SPECIFIC GRAVITY: (20 °C) N/A
SOLUBILITY IN WATER: (20 °C) Soluble

SECTION X: STABILITY AND REACTIVITY

SUBSTANCES TO BE AVOIDED: N/A
HAZARDOUS, COMBUSTION, OR DECOMPOSITION PRODUCTS: N/A

SECTION XI: TOXICOLOGICAL INFORMATION

TOXICITY DATA: N/A
INHALATION: May be harmful by inhalation.
EYE CONTACT: May cause eye irritation.
SKIN CONTACT: May cause skin irritation, may be harmful by skin absorption.
INGESTION: May be harmful if swallowed.
PROLONGED EXPOSURE: N/A
CHRONIC EFFECTS: N/A
RTECS NUMBER: N/A
ADDITIONAL INFORMATION: THE CHEMICAL, PHYSICAL, AND TOXICOLOGICAL PROPERTIES HAVE NOT BEEN THOROUGHLY INVESTIGATED. Additional harmful properties cannot be ruled out. The product should be handled with the normal caution accorded chemicals.

SECTION XII: ECOLOGICAL INFORMATION

N/A

SECTION XIII: DISPOSAL CONSIDERATIONS

There are no uniform regulations for the disposal of chemicals or residues. Dispose of container and unused contents in accordance with federal, state and local requirements.

SECTION XIV: TRANSPORT INFORMATION

DOT: None of the components are regulated.

SECTION XV: REGULATORY INFORMATION

SARA: None of the components are regulated.

SECTION XVI: OTHER INFORMATION

DATE OF PREPARATION: March 1, 2015.
DISCLAIMER: For research use only. The above information is believe to be correct but does not purport to be all inclusive and should be used only as a guide. MultiTarget Pharmaceuticals shall not be held liable for any damages or other consequences resulting from handling or from contact with the above product.